


ALERGENI NA LICU MJESTA

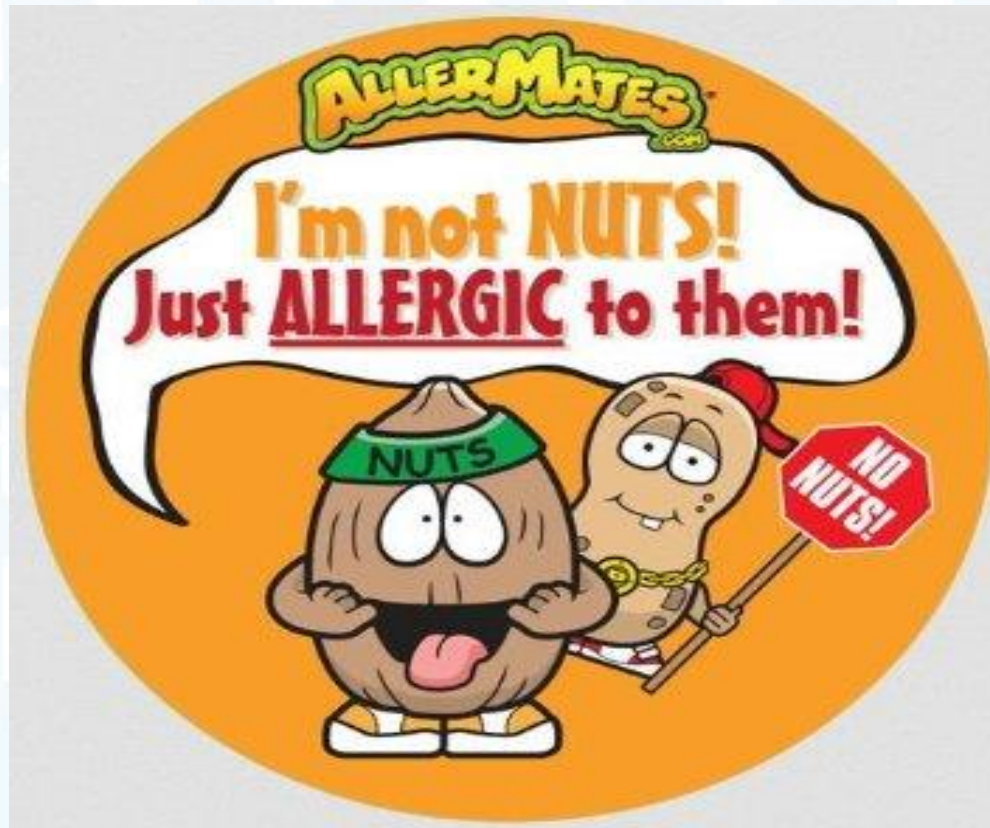

Koprivnica, 29.09.2016.

REVEAL MULTI TREENUT



...a Vi ?

REVEAL MULTI TREENUT



**Jedinstveni laterl flow test
detektira pojedinačni ili
kombinaciju od šest orašastih
plodova:**

Badem

Lješnjak

Pekan Orah

Indij.oraš.

Pistachio

Vrijeme testiranja: 10 min

Broj testova: 10 uzoraka/briseva

**Resultati: 2 testne
linije**

**1 linija negativno
2 linija pozitivno**



Reveal Multi-Treenut Postupak

Ostaci ispirnih otopina- ekstrakcija



1. Otvorite vrećicu sa puferom i dodajte u uprjetu za uzorak.



2. Dodajte 1 mL uzorka.




3. Zatvorite i protresite 1 min.

Reveal Multi-Treenut Postupak

Brisevi okoliša-ekstrakcija



1. Otvorite vrećicu sa puferom i dodajte u epruvetu za uzorak

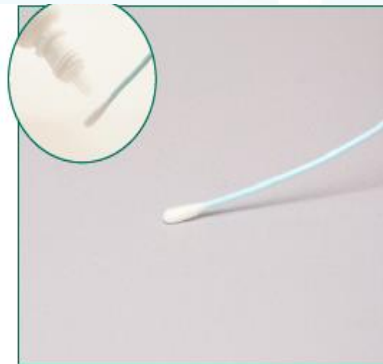

2. Uzmite uzorak-bris sa površine.



3. Vratite bris u eprivetu sa puferom i prelomite ga.



4. Zatvorite i protresite 1 min.

Reveal Multi-Treenut postupak

Testiranje uzorka



1. Otvorite epruvetu sa uzorkom i napunite čep sa ekstraktom iz epruvete.

2. Umočite Reveal test u čep tako da se otvor testa ipuni tekućinom.

3. Ostavite tako dugo dok vidite da tekućina prolazi kroz testni prozorčić.

4. Stavite test na ravnu površinu i ostavite 10 minuta.

Reveal Multi-Treenut Rezultati



← **NEGATIVNO**

← **POZITIVNO**

VAŽNO: linija je linija; pojava linije bilo kakvog intenziteta na T i C poziciji indicira POZITIVAN rezultat

Reveal Multi-Treenut

Brisevi okoliša 5 - 40 $\mu\text{g}/100\text{cm}^2$

Treenut	Sensitivity ($\mu\text{g}/100\text{cm}^2$)		
	Stainless Steel	Teflon	Plastic
Walnut	5	5	5
Hazelnut	20	10	20
Cashew	20	20	20
Almond	20	20	20
Pistachio	40	20	40
Pecan	20	10	10

CIP otopina – 5-10 ppm

CIP otopina LOD 5-10 ppm za svaki od navedenih orašastih plodova .

Napomena: prisustvo detergenta u ispiranom ostatku može utjecati na LOD.

HVALA !

